

## SHOC INSTRUMENT CHANGE NOTES: APPLICABLE TO 40" TELESCOPE ON

May 2012	Jaci Cloete
March 2014	Willie Koorts
July 2014	Jaci Cloete
Feb 2015	Piet/Jeran
May 2015	DBC
Nov 2015	Keegan Titus
Dec 2015	Michael Rust
Jan 2016	Willie Koorts
Feb 2016	Willie Koorts
6 Sept 2016	Dave Carter
19 Oct 2016	Carlson Mulaudzi
01 Mar 2018	Avhaphani Mulaudzi
08 Oct 2018	Avhaphani Mulaudzi

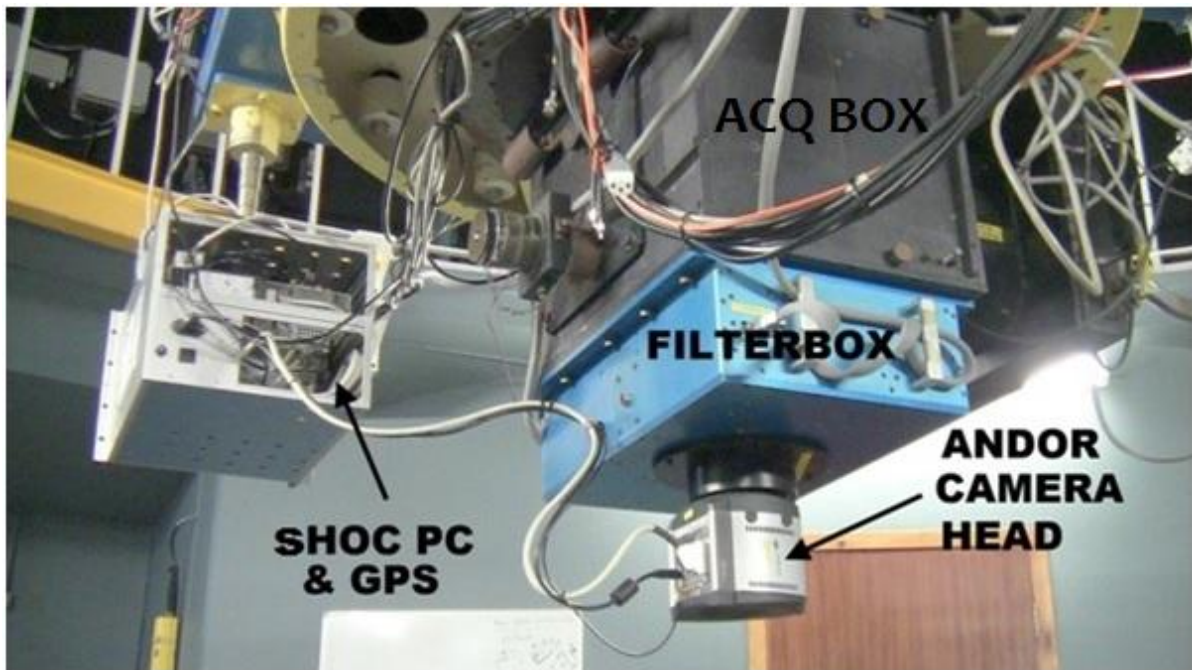


Figure 1a: SHOC on the 40". View from the NW

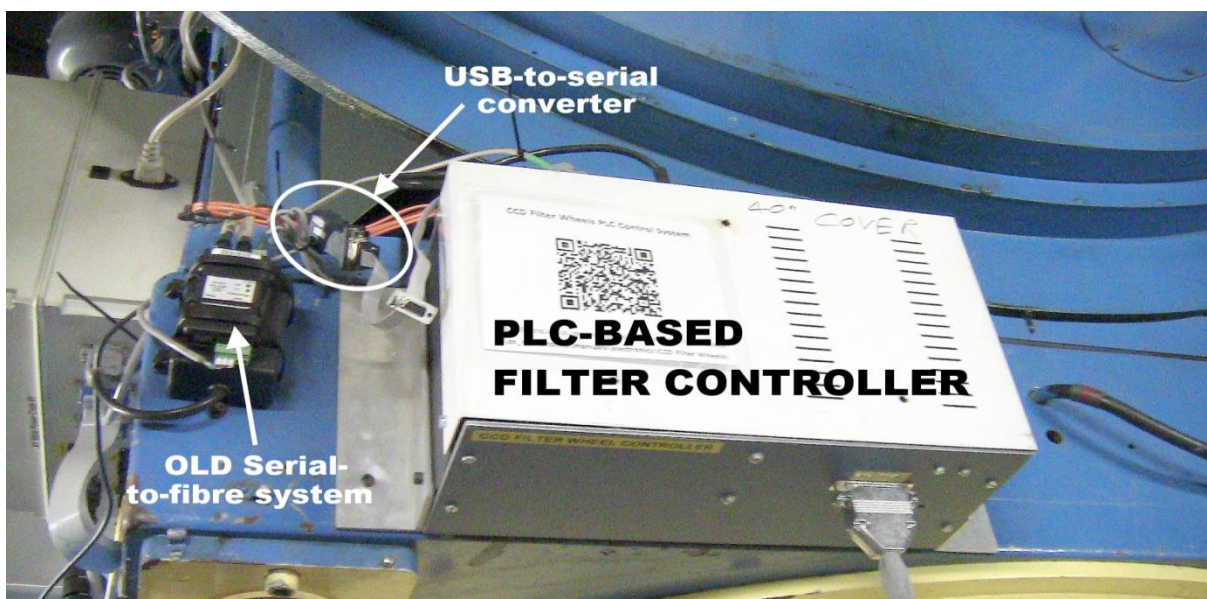


Figure 1b: Filter wheel controller

### In the Warmroom:

1. Ensure that **TCS** is locked out by turning the **TCS Lockout** switch on the mimic panel (refer to figure 2) to **ON** state. TCS Control indicator will flash as a warning that the TCS is locked out.



**Figure 2: TCS LOCKOUT SWITCH**

### ON THE TELESCOPE:

1. Mount SHOC PC Crate on Telescope (see picture). See Appendix for the implications of mounting the "wrong" PC Crate. Do NOT Connect the mains power cable at this stage.
2. Ensure that XY slides and autoguider are mounted on the acquisition box (permanently mounted on telescope) on the South side.
3. Ensure that IEC mains cable and 15-way ribbon cable are connected to the acquisition box.
4. Mount Andor Camera Unit on Telescope – see picture. (Ensure Camera lens has been cleaned by gently blowing with compressed air.)
5. Ensure Check that the correct filter wheels are installed in the filter box – refer to the 'Filter Rota' list on the electronics web page.
6. Connect the Filter Wheel Controller (permanently mounted on telescope) filter cables to filter wheels. Connect the USB-to-serial cable from filter wheel controller to SHOC PC (see pic.)
7. Connect GPS Antenna (permanently mounted on telescope) to GPS Unit inside PC Crate
8. Connect Network Cable (transceiver permanently mounted on telescope) to PC Crate
9. Connect SHOC PC to Andor Camera Unit
10. Connect PSU to Andor Camera Unit

11. Connect GPS trigger signal cable to Andor Camera unit “ext trig”
12. Connect PC Crate to Telescope mains (normal Kettle plug lead)
13. Switch on Crate Power & SHOC PC (in front behind black tape)

#### IN THE WARMROOM:

1. Useful info may be available at: <http://topswiki.saao.ac.za/index.php/SHOC> - especially if there’s any doubt as to whether these instrument change notes are up-to-date with possible software updates . . .
2. Run the appropriate script as follows:
  - 3.1 Open xterminal, type in “ssh [ccd@shoc40in.suth.saao.ac.za](http://ccd@shoc40in.suth.saao.ac.za)” when prompt for a password type “**Saaoccd**”.
  - 3.2 Run **shocboxswitch.sh** command, then when asked “Do you wish to revert to the default?” type “**NO**”
  - 3.3 When asked “**which telescope should be used?**” then type in **40in** and when prompt for a password type “**Saaoccd**” the script will run.
3. Connect to the web browser: <http://shoc40in.suth.saao.ac.za:5000>. When prompted to log in, use the username and password below:
 


**\*VERY IMPORTANT!** See Appendix for a detailed explanation by Amanda about the implications of mounting the “wrong” SHOC Computer on the “wrong” Telescope.

Username: shoc40in

Password: Saa040in

There are three tabs, for **Filter control**, **GPS control**, and **Camera control**

4. Go to the **Camera** tab, click the “**Turn Camera on**” button – see figure 3. If the camera switches on OK the “Control” tab GUI will appear – See Figure 4.
- 5.

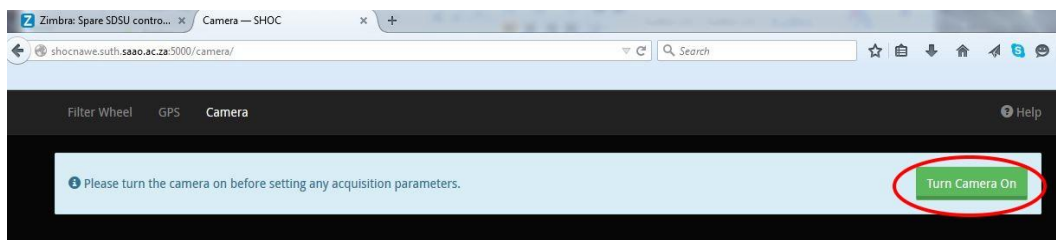

**Figure 3: Initial state of "Camera" tab**

6.


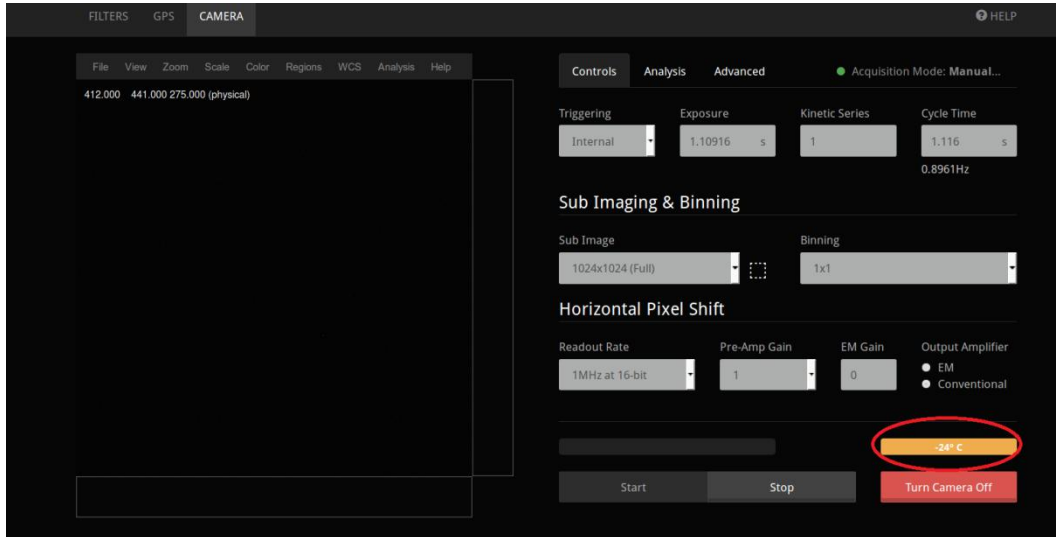


Figure 4: Camera "Control" tab GUI

m

era temperature: See the temperature display/button on the bottom RH corner. It should have orange background and the reading should start dropping.


- a. Click on the Temperature display button. The "advanced" tab GUI should appear – see Figure 5.
- b. To avoid unnecessary stress to the thermoelectric cooler, click on the "Set Temperature" box (Figure 5) and change the set point to -25°C. Check that the temperature settles at the new set point. Note that the servo seems to overshoot quite a lot. Once at the set point, the button background colour changes to green. This can take a few minutes . . .

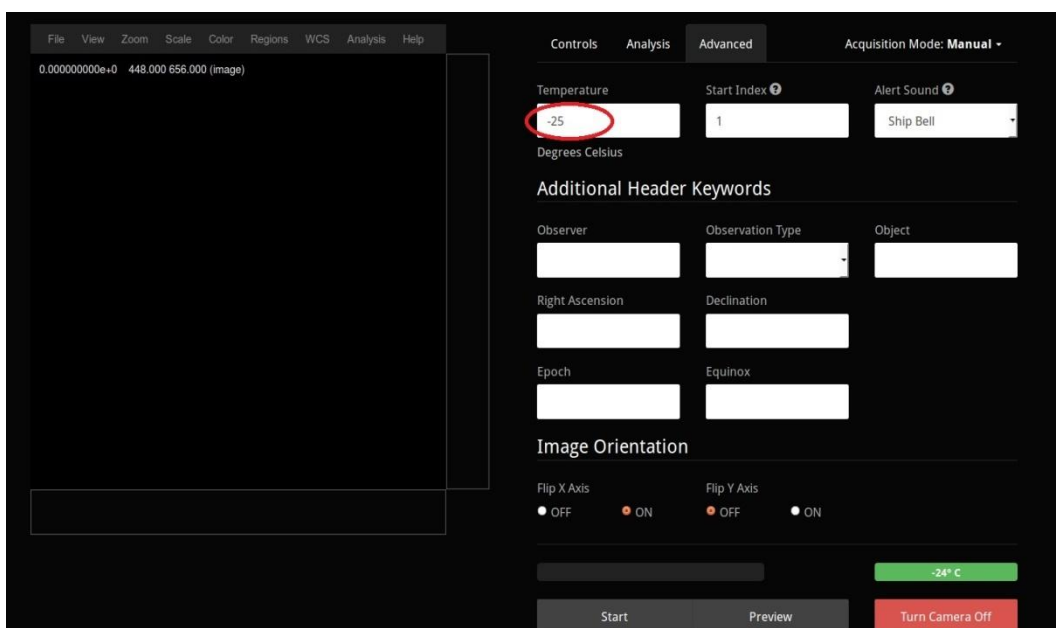

**Figure 5: "Advanced" tab GUI**

7. 

Check that the camera is reading out: Click on the "Controls" tab (Figure 6), confirm the setting is full frame, conventional mode (the default) and press "preview". Images should appear every ~ 1.2 sec. Once confirmed that images are coming through, press "stop". Note that camera has no shutter, therefore before bias frames can be taken, the dome lights must be switched off, ensure dome shutters and cover is closed (it's not necessary that it should be completely dark inside the dome, but there should be no direct light falling on the optical path). Running the cursor over the image display area will bring up a notification panel with X,Y coordinates and the count at this position. The counts should be around 400 or so.

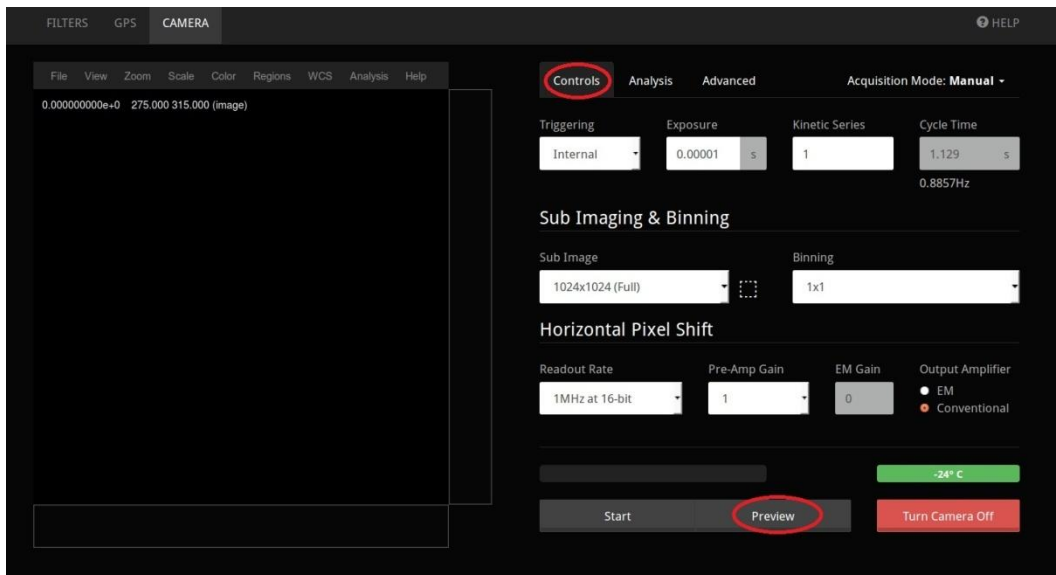


**Figure 6: Camera "Controls" Tab**

8. Check that the camera fan is running. (Check this at the camera)
9. Turn the camera off: Click on the "Turn Camera Off" button at the RHS bottom corner of the GUI.


Check the Filter & GPS operation: Click on the "Filter Wheel" and "GPS" Tabs at the top RH corner of the web GUI to access the controls. See Figures 7 & 8.

10. On the web page Filter Tab initialize and move filters to check their operation.

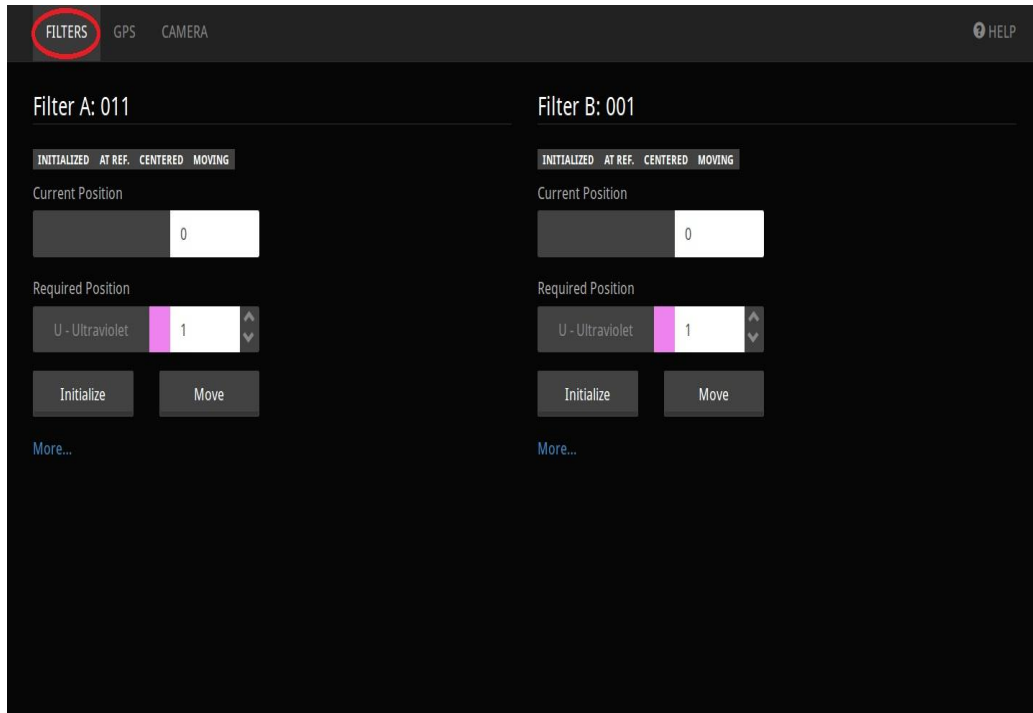

**Figure 7: Filter Wheel GUI**

11. On the web page GPS Tab, just check to ensure that there is no Antenna Fault (If this happens check that the antenna has been connected properly to the GPS unit in the port labelled "antenna" also refer to page 18 of the user manual under the section titled "hardware fault monitoring" point 2.)

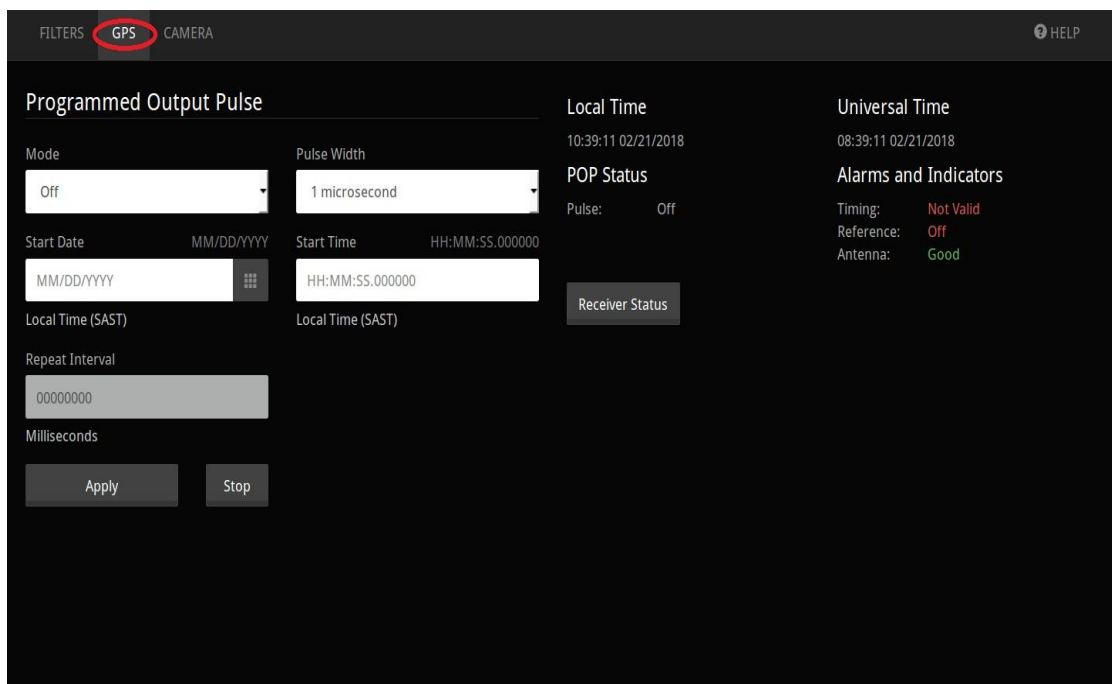

**Figure 8: GPS Tab GUI**



12. GPS data acquisition is an automatic process and will start as soon as dome is opened.
13. Do the usual checks on the TCS/XYslide/acquisition camera hardware.
14. **NB!!! Ensure that TCS Lockout switch on the mimic panel is set to ON state.**

Further information and user manuals (for trouble shooting) can be found at <http://shoc.sao.ac.za/Documentation.html>

---

### **Appendix:**

In an email dated 15 January 2016, Amanda explained the implications of mounting the “wrong” CHOC Computer on the “wrong” telescope, as follows:

“If the boxes are switched (i.e. shocnawe goes on the 40in or shocndisbelief goes on the 74in), then the observer will not be able to connect to the machine using the appropriate website and account (which is shoc40in or shoc74in for each telescope, respectively).

The instrument could be accessed by logging into the WRONG account, so if shocndisbelief were mounted on the 74in, then the 74" observer could run the shoc40in web interface and transfer data as shoc40in. This would be functional, but then all of the data paths would be incorrect because files would be stored under the 40in telescope folders, and the observer is likely to get confused when logging in and out. Things could get really hairy in this case if for some reason multiple SHOCs were mounted and an observer mistakenly logged into the wrong one, thus controlling the instrument on a different telescope.

The preferred solution is for IT to know how to change whatever information is required so that the SHOC box can be correctly tied to the correct telescope username where it is mounted.”